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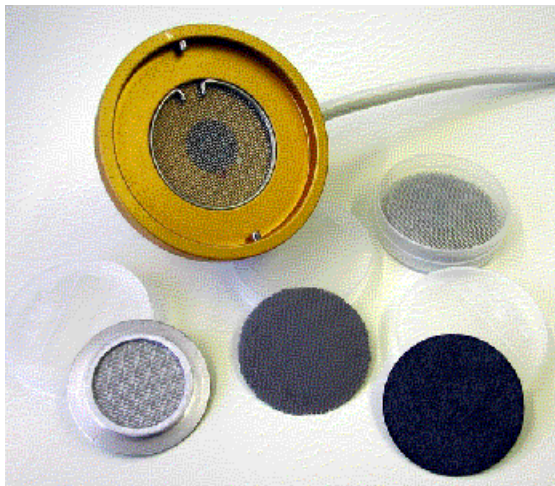
These specially designed filters shall be used to provide protection against chemical perturbation. Mineral acids, halogens, powerful oxidizing agents and corrosive substances may damage the cells and reduce their lifetime. Chemical filters shall be replaced, when a recalibration of the cell becomes necessary more often, when replacing a cell, but latest after one year. The eVALC needs not to be replaced, if not mechanically damaged.

The higher the concentration of an aggressive substance, the less effective will be the protection capability of a chemical filter, so the cell lifetime still can be reduced. If the concentration exceeds certain limits, no protection is guaranteed.

-> Only practical testing can answer the question about exchange cycles of a filter.

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Novasina FILTER SYSTEMS

eVC-21 Chemical Filter

acetic acid (eg. mustard)
formic acid
butyric acid
lactic acid
sulphuric acid
sulfur dioxide
hydrogen sulfide
chlorine
carboxylic acids, aliphatic
hydrogen halides
hydrogen peroxide
formaldehyde
jodine
chlorine, hypochloric acid

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eVC-26 Universal Filter

oil fog
fine dusts
amines
ozone
nicotine
aromatic hydrocarbons
perchlor-, trichlor-ethylene
chloroform
nitrogen oxides

Art.nr. 1111003

Oelnebel
Feinstaub
Amine
Ozon (O₃)
Nikotin
arom. Kohlenwasserstoffe (Benzol, Kerosen, Toluol)
Perchlor-, Trichlorethylen (C₂HCl₃)
Chloroform (CHCl₃)
Stickoxyde (NO_x)

Redox Filter

primary, secondary alcohols
glycerol
ethanol (ethyl alcohol)
isopropanol (isopropyl alcohol)
propanol (propyl alcohol)
propylen-glycol
polyethylen-glycol
methanol
phenol

Art.nr. 1117212

primäre, sekundäre Alkohole (RCH₂OH, R₂CHOH)
Glycerin [CH₂(-OH)-CHOH- CH₂(-OH)]
Ethanol (C₂H₅OH)
Isopropanol [(CH₃)₂CHOH]
Propanol (C₃H₇OH)
Propylen-Glykol (CH₃CHOH CH₂OH)
Poly-Ethylen-Glycol (PEG)
Methanol (CH₃OH)
Phenol (C₆H₅OH)

yeast and hop aromas
coffee aromas
artificial fruit aromas
onions, garlic, horseraddish
furanes (bread aromas)
ketones (eg. acetone or MEK)
partly *amines*
perfumes, flavours, fragrances

Hefe-, Hopfenaromen
Kaffeearomen
künstliche Fruchtaromen
Zwiebeln, Knoblauch, Meerrettich
Furane (Brotaromen)
Ketone (zB. Aceton od. MEK) {(CH₃)₂CO, (CH₃)CO(C₂H₅)}
teilweise *Amine* (NH₂R, NHR₂, NR₃)
Parfüms, Duftstoffe

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